

## How to install the BarTender software

Please skip directly to step 3 if you purchased a full license of the BarTender software

1. Please go to:

<https://www.seagullscientific.com/30-day-trial/>

and sign up for a free user portal account.

2. Once logged in, please navigate to **Downloads** to get the license code for your 30-day trial.

3. Locate the software installation file on this flash drive and start the installation process.

4. Follow the instructions on screen to install the software.

5. When prompted, please enter the license key code into the software in order to activate it. (Please note that you will need an internet connection to do so).

## How to install Cody-3 driver

To install the required printer driver, locate the “**Generic\_2021.2\_M-2**” file in the “Drivers” folder to start the installation process.

Please contact us for technical support at: [support@ga-international.com](mailto:support@ga-international.com)

or call us toll free at **1-800-518-0364 (US and Canada).**

**+1 (450) 473-9420 (International).**

**ga-international.com**

**Labtag.com**